

## OPERATING INSTRUCTIONS INDUCTFLOW CHAMBERS



1. Place the chamber on a flat surface with the ports positioned away from the operator. Verify floor liner is in place.
2. Attach the Fresh Gas adapter on the bottom (red) port and the Waste Gas Adapter on the top (blue/black) port.
3. Turn Waste Gas Evacuation system ON and/or verify W.A.G. filter is connected.
4. Anesthetize the animal following standard protocol.
5. Once the animal is fully anesthetized, turn the vaporizer dial off and allow several seconds for the oxygen/carrier gas flow to aid in reduction of anesthetic vapor left within the chamber (passive style chamber only).
6. Gently tilt the chamber to safely position the animal to the far opposite side (port side) of the chamber.
7. While keeping the chamber facing away from the operator, slide the lid open only far enough to allow safe removal of the animal. Immediately close the lid to prevent escape of remaining waste anesthetic gas.
8. When use of induction chamber is complete, continue flowing oxygen/carrier gas through the chamber and/or keep evacuation system on to allow all remaining anesthetic to be removed.

### **CLEANING:**

Clean with mild soap and water following standard protocol.

**SOMNI PROVIDES UNPARALLELED CUSTOMER SERVICE, CLINICAL AND TECHNICAL SUPPORT.**

For more information or pricing, please contact our sales team.

☎ 1-877-637-3625    ✉ info@somniscientific.com

🌐 www.somniscientific.com



*Proudly machined, assembled  
and calibrated in the USA*